BOD₅ Glucose – glutamic acid standard

Principle

LCA 555 *addista* is a control standard system for cuvette test LCK 555 BOD₅. The standard contains glucose and L-glutamic acid.

Storage Information

The test reagents are stable at +2 to +8°C up to the expiry date given on the package.

Special note

Always prepare fresh standard and discard after use. Carry out the analysis in duplicate if possible.


Characteristic Data

Recovery rate for optimal inoculation	205 +/- 45 mg/l BOD ₅
---------------------------------------	----------------------------------

Safety Advice

On grounds of quality and reliability, the analysis should be carried out only with original Dr. Lange accessories.

Cuvettes with freeze-dried contents contains: Natriumnitrat
R: 8-22-36 S: 22-24-41

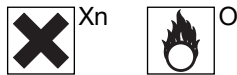R: (8) Contact with combustible material may cause fire. (22) Harmful if swallowed. (36) Irritating to eyes.

S: (22) Do not breathe dust. (24) Avoid contact with skin. (41) In case of fire and /or explosion do not breath fume.

Procedure/Evaluation **LCA 555**

BSB₅ Glucose - glutamic acid standard

Edition 98/04

1. Prepare the standard solution

Into a cuvette containing freeze-dried contents, pipette
Distilled water 5.0 ml

Close the cuvette and shake vigorously. After **3 min pipette 2.2 ml of this solution** into a 50 ml measuring flask, top up with distilled water and shake vigorously.

2. Analysing the standard

Standard solution (50 ml measuring flask) 0.5 ml
directly into sample cuvette LCK 555 pipette.
Top up the sample cuvette with dilution water.

3. Evaluation

After **5 days, evaluate in measuring range B2** (50 - 275 mg/l) against the dilution water cuvette that was prepared in parallel.

Note

The theoretical value of 205 +/- 45 mg/l corresponds to the generally accepted scatter of the biological measuring procedure for BOD₅.

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