T1
2000 mg/L: K ⁺
1500 mg/L: Na ⁺
1000 mg/L: Cl ⁻
500 mg/L: CZV / COD *)
250 mg/L: Ca ²⁺
100 mg/L: Ag ⁺
50 mg/L: Pb ²⁺ , Zn ²⁺ , Ni ²⁺ , Fe ³⁺ , Cd ²⁺ , Cu ²⁺
20 mg/L: Fe ²⁺
10 mg/L: Co ²⁺
5 mg/L: Cr ⁶⁺
2 mg/L: NO ₂ ⁻

*) (Kaliumwaterstofftalaat)
(Potassium hydrogen phthalate)

Datatablel / Data table

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GB LCK 340 Nitrate

Please check the "Edition Date" (see data table) and read the "Note". Safety advice and expiry date on package.

Principle

Nitrate ions in solutions containing sulphuric and phosphoric acids react with 2.6-dimethylphenol to form 4-nitro-2.6-dimethylphenol.

Range of Application

Waste water (beware of interferences!), drinking water, raw water, surface water, soils, substrates, nutrient solutions

Interferences

The ions listed in **T1** have been individually checked up to the given concentrations and do not cause interference. We have not determined cumulative effects and the influence of other ions. High loads of oxidizable organic substances (COD) cause the reagent to change colour and give high-bias results. The test can thus only be used for waste water analyses if the COD is less than 500 mg/L.

The measurement results must be subjected to plausibility checks (dilute and/or spike the sample).

Removal of Interferences

Nitrite concentrations of more than 2.0 mg/L interfere (high-bias results) and can be removed by the addition of a spatula-tipfull of amidosulphonic acid. The chloride can be precipitated out as silver chloride by adding silver sulphate.

High calcium concentrations cause turbidity. This interferes with the determination but can be prevented by adding a spatula-tipfull of EDTA to the sample.

pH sample3 – 10
Temperature sample/reagents20 – 24°C

In case of not working at the right recommended temperature an incorrect result may be obtained.

Not more than 3 hours should elapse between sampling and analysis. **Store in a cool place!**

Note

Change of factor for all types of photometers (except LASA aqua).

T1
2000 mg/L: K ⁺
1500 mg/L: Na ⁺
1000 mg/L: Cl ⁻
500 mg/L: CSB / DCO / COD *)
250 mg/L: Ca ²⁺
100 mg/L: Ag ⁺
50 mg/L: Pb ²⁺ , Zn ²⁺ , Ni ²⁺ , Fe ³⁺ , Cd ²⁺ , Cu ²⁺
20 mg/L: Fe ²⁺
10 mg/L: Co ²⁺
5 mg/L: Cr ⁶⁺
2 mg/L: NO ₂ ⁻

*) (Kaliumhydrogenphthalat)
(Hydrogènthalate de potassium)
(Potassio idrogenoftalato)

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